


# ACID RESIN

MILESTONE APPLICATION NOTE FOR ACID DIGESTION  
SK-CHEMICALS-004

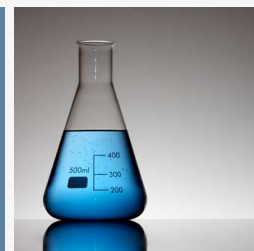

## SUMMARY

This method represents a guideline for the microwave acid digestion of Acid resin samples with the Milestone ETHOS UP and ETHOS EASY systems. This procedure may need to be optimized to match your specific analytical goals.

## INSTRUMENTATION

Milestone ETHOS UP or ETHOS EASY microwave units, SK-15 high-pressure rotor, T2 easyTEMP temperature control. Options T1 temperature control and P1 or P2 pressure control.

## SAMPLE WEIGHT AND REAGENTS

|               |                                                             |
|---------------|-------------------------------------------------------------|
| Sample weight | Up to 500 mg                                                |
| Reagents      | 9 mL HNO <sub>3</sub><br>1 mL H <sub>2</sub> O <sub>2</sub> |



## MICROWAVE PROGRAM

| Step | Time     | T2    | Power*  |
|------|----------|-------|---------|
| 1    | 00:20:00 | 210°C | 1800 W* |
| 2    | 00:15:00 | 210°C | 1800 W* |

### \* POWER SETTINGS

The power of the system must be set according to the table below:

|                   |        |
|-------------------|--------|
| 3 or less vessels | 800 W  |
| 4-8 vessels       | 1200 W |
| 9-15 vessels      | 1800 W |

## GENERAL PRECAUTIONS

Always use hand, eye and body protection when operating with the microwave system.

Subject to change without notice.  
For additional information  
please visit [www.milestonesrl.com](http://www.milestonesrl.com)  
or contact [application@milestonesrl.com](mailto:application@milestonesrl.com)

